

Using the Macintosh Data Recorders with Logger Pro

Setting up the Data Recorder

1. Remove the data recorder and any accessory materials from the storage location. The Vernier LabPro box and AC power supply adapter may be located with the data recorder or with the Vernier supplies for the laboratory.
2. Check that all connections are secure:
 - a. Be sure the power cable is connected to the data collector, and that the power cable is plugged into a power supply
 - b. Connect the cable between USB port on the data collector and the Vernier Labpro box. Be sure the LabPro is also connected to its power supply.
3. Plug in the sensor to be used into the Labpro "Ch1" (on the side of the LabPro).
4. Start the data collector.


Interface Cable

Most of the LoggerPro peripheral devices (sensors) need to have an interface cable to connect to the Vernier LabPro box. If one is not provided with the laboratory supplies or with the data recorder check with your instructor.

Setting Data to Read for 5 Decimal Places for Labs Using the Gas Pressure Sensor

1. Choose Data Menu
2. Choose Column Options
3. Choose Pressure
4. Set to 5 decimal places

Using the Data Recorder's Trackpad

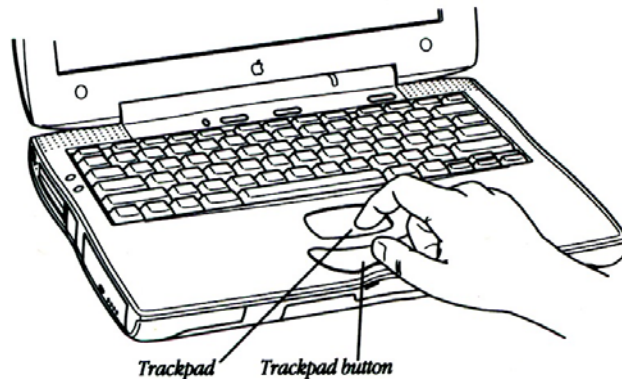

To navigate with the data recorder's trackpad, hold down the trackpad button with your thumb and move your finger on the trackpad to "drag" the pointer on the screen to the desired location. Release the trackpad button. The cursor should be in the location you want.